
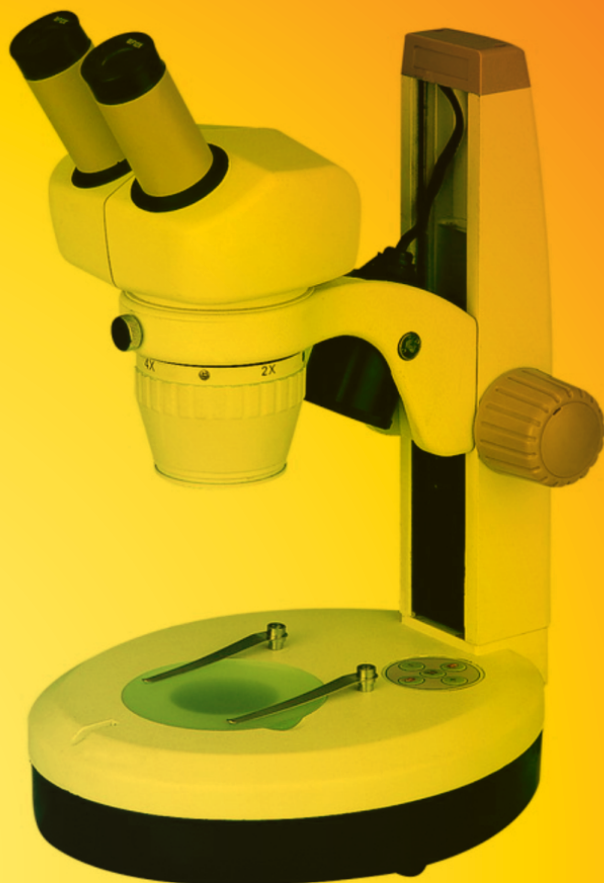
# Levenhuk 5ST stereo microscope

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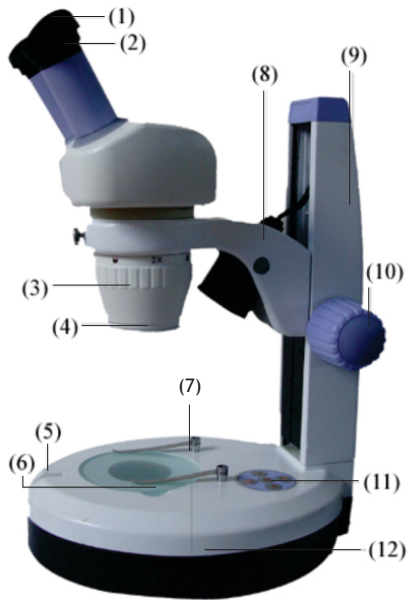**levenhuk**   
Zoom&Joy

- 1 Rubber eyecup
- 2 Eyepiece
- 3 Revolving nosepiece
- 4 Secondary objective lens adapter
- 5 Locking screw
- 6 Stage
- 7 Stage clips
- 8 Upper illumination source
- 9 Pole
- 10 Focus knob
- 11 Illumination switch
- 12 Base

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## Specifications and the kit

Eyepiece	10x
Objective lens	2x, 4x
Interpupillary distance range	55 mm - 75 mm (2.17-2.95 in)
Microscope head inclination angle	60°

Magnification chart:

Secondary objective lens	Objective lens		1x	2x	2x	2x
	Eyeiece					
0.5x	10x		5x	10x	15x	20x
2x			20x	40x	60x	80x
0.5x	15x		7.5x	15x	22.5x	30x
2x			30x	60x	90x	120x

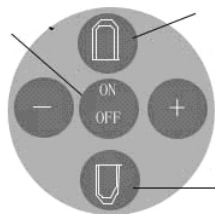
The following accessories are not included in the kit and have to be purchased separately: 15x eyepiece; 1x, 3x objective lenses; 0.5x, 2x secondary objective lenses

## Using the microscope

**Illumination.** You can turn on both illumination sources by pressing the ON/OFF button. By pressing the '+' and '-' buttons you can set the brightness level.

**Note:** you can turn on either the upper or the lower illumination source individually by pressing the corresponding buttons.

Illumination power button

Upper illumination source switch

Lower illumination source switch

- Place a slide on the stage and lock it in place with stage clips.
- Adjust the interpupillary distance by holding both eyepieces with your hands and bringing them closer or moving them apart until a double image turns into a single image.
- Perform diopter adjustment, if necessary, and focus your view.
  - (1) Look through the right eyepiece. Focus the view by rotating the focus knob.
  - (2) Look through the left eyepiece. Focus the view by rotating the diopter adjustment ring.

## Caution

**Never, under any circumstance, look directly at the Sun, another bright source of light or at a laser through this device, as this may cause PERMANENT RETINAL DAMAGE and may lead to BLINDNESS.**

Take necessary precautions when using the device with children or people, who have not read and fully understood these instructions.

Do not try to take the device apart on your own. For repairs of any kind, please contact your local specialized service center.

## Care and maintenance

- After unpacking your microscope and before using it for the first time check for integrity and durability of every component and connection.
- Do not touch the optical surfaces with your fingers. You can use a soft napkin or a cleaning wipe, dipped in ether or absolute alcohol, to clean lenses. We recommend using special optics cleaning fluids from Levenhuk. Do not use any corrosive or acetone-based fluids to clean the optics.
- Abrasive particles, such as sand, should not be wiped off lenses, but instead blown off or brushed away with a soft brush.
- Do not apply excessive pressure when adjusting focus. Do not overtighten the locking screws. Protect the device from sudden impacts and excessive mechanical force.
- Store the device in a dry, cool place away from dust, hazardous acids and other chemicals, away from heaters, open fire and other sources of high temperatures. Do not use the device for lengthy periods of time, or leave it unattended in direct sunlight. Keep the device away from water and high humidity.
- Be careful during your observations, always replace the dust cover after you are finished with observations to protect it from dust and stains.
- If you are not using your microscope for extended periods of time, store the objective lenses and eyepieces separately from the microscope.

## International Warranty

- The Vendor guarantees that the quality of the Levenhuk product purchased complies with the technical documentation requirements on conditions that the consumer observes rules of transportation, storage conditions and operating instructions.
- Levenhuk Ltd. warrants the product against defects in materials.
- During the warranty period the Buyer can return the defect product to the Vendor or to Levenhuk service center. Levenhuk Ltd. will repair or replace the defect product at its own discretion.
- No claims are accepted in case the properly filled warranty slip is absent or contains corrections, or if the defect product has not been provided by the Buyer. The manufacturer or the seller is not liable for any damages caused by misuse of the product. No guarantee is provided in cases the product was used for purposes other than that intended or has mechanical damages, scratches, cracks, optics damages, as well as in cases the product is broken-down due to impact damages, squeezing or stretching or the product has been disassembled or repaired by unauthorized personnel.
- This warranty does not cover components and/or accessories with limited use period, batteries, etc.

The period of warranty is 3 (three) years beginning on the date of purchase. Please keep the warranty slip along with your receipt.

For any warranty services please contact the local Levenhuk branch.

**Levenhuk Worldwide:** USA: [www.levenhuk.com](http://www.levenhuk.com), Czech Republic: [www.levenhuk.cz](http://www.levenhuk.cz), Russia: [www.levenhuk.ru](http://www.levenhuk.ru), Ukraine: [www.levenhuk.com.ua](http://www.levenhuk.com.ua), EU: [www.levenhuk.eu](http://www.levenhuk.eu)

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